


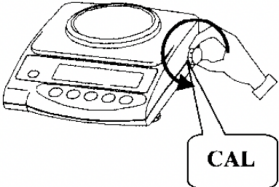
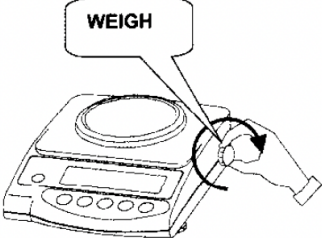


Instruccions calibratge balança KERN EW/EG-N

Adjustment procedure:
Check that the surrounding conditions are stable.
A short warm-up time of about 10 minutes is recommended for stabilisation.

Operation	Display
Switch-on balance with button  , after a short time [S.A. CAL] will appear.	<div>S.A. CAL</div>
Press buttons  and  at the same time and release at the same time, [Wait] is briefly displayed.	<div>Wait</div>
After that [CAL.0] will appear flashing, Zero point will be saved.	<div>CAL. 0</div>
After that [CAL.on] will appear.	<div>CAL. on</div>
Rotate the turn knob on the right side of the balance in to position CAL.	<div>CAL. on</div>
 Adjustment is carried out automatically. [CAL.on] will appear flashing.	
The display changes automatically from [CAL.on] to [CAL.off] The process of adjustment is completed.	<div>CAL. oFF</div>
Turn the adjusting knob on the right side of the balance to the WEIGH position.	<div>buSY</div>
 Calibration is now finalised. The balance will automatically return to weighing mode.	<div>End.</div>

