

DRYING SOLID SAMPLES WITH THE VACUUM LINE

Alicia Mate González & Toni Durán Mota

MATERIAL AND PREPARATION OF THE ASSEMBLY

put the sample
inside the
schlenk for
drying samples



vial with sample to be dried

cover the vial
with parafilm
and make a few
holes with the
needle



close the
system with
the screw cap



*check that the
screw caps have
their adapter
well placed
before closing
the system*

CONNECT THE ASSEMBLY TO THE VACUUM LINE
WAIT UNTIL THE SAMPLE IS DRIED (15-20 MINUTES)

